


Infratec™
Installation Guide
6008 9468 / Rev. 1


1

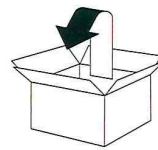

x1



x1

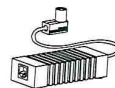

x1




x1




x1





x1

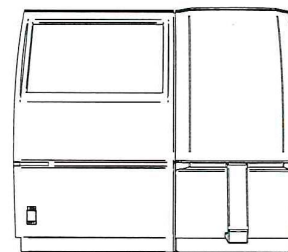

4

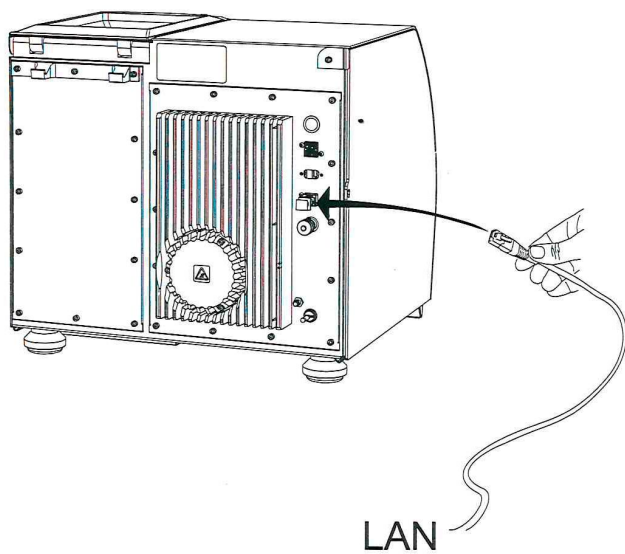


5

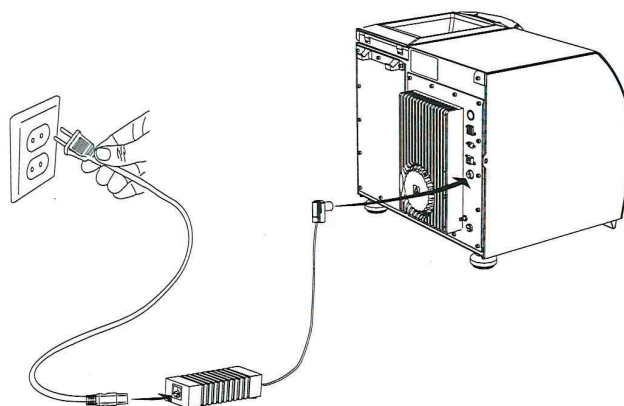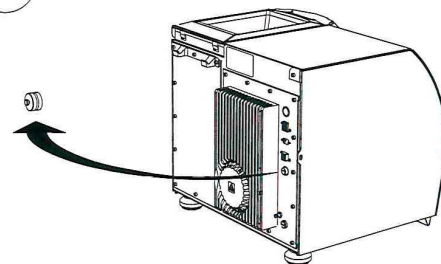


100-240 VAC




2

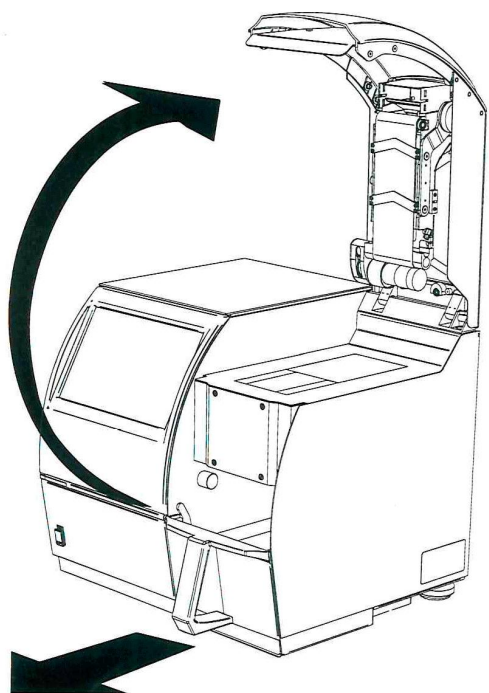

3

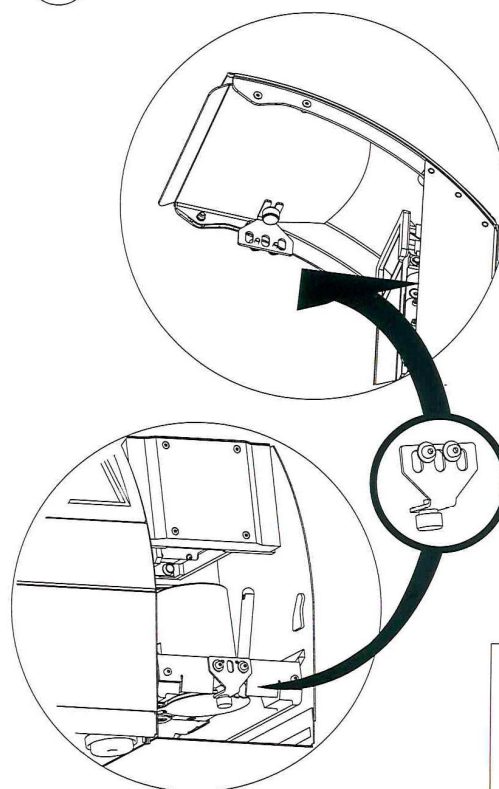

6


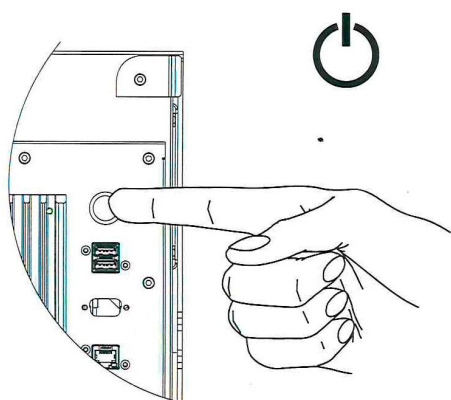


7



-  Installation Guide
-  Owners Guide
-  Quick Guide
-  Safety Manual
-  Spare Parts Manual
-  User Manual

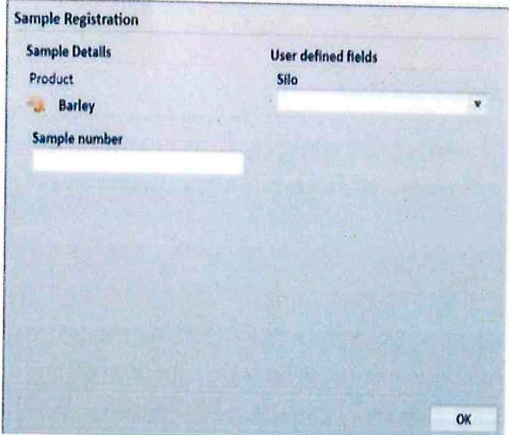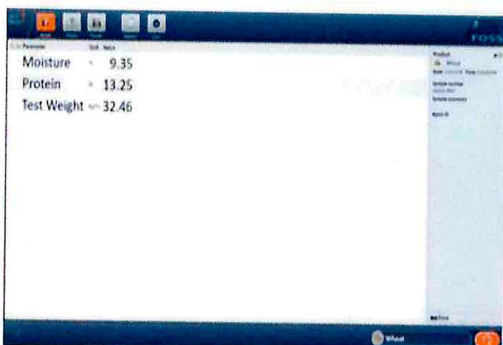

3. Press Start button.	
4. Register sample number and other relevant sample information after start of analysis.	
5. Once the analysis is complete, the result will be presented on the screen and the sample is collected in the drawer.	

Turning Off the analyzer

1. Go to the Care Menu.
2. Press SHUT DOWN.
3. Press OK.

Quick Guide

Infratec™ **6008 9464 / Rev. 1**



Please read safety precautions and warranty statement in the Safety Manual prior to use.

Before operating the instrument, ensure that the instrument and all external equipment are installed and connected according to the User Manual.

Note that your installation may differ from the one described.

ISIsScan Nova

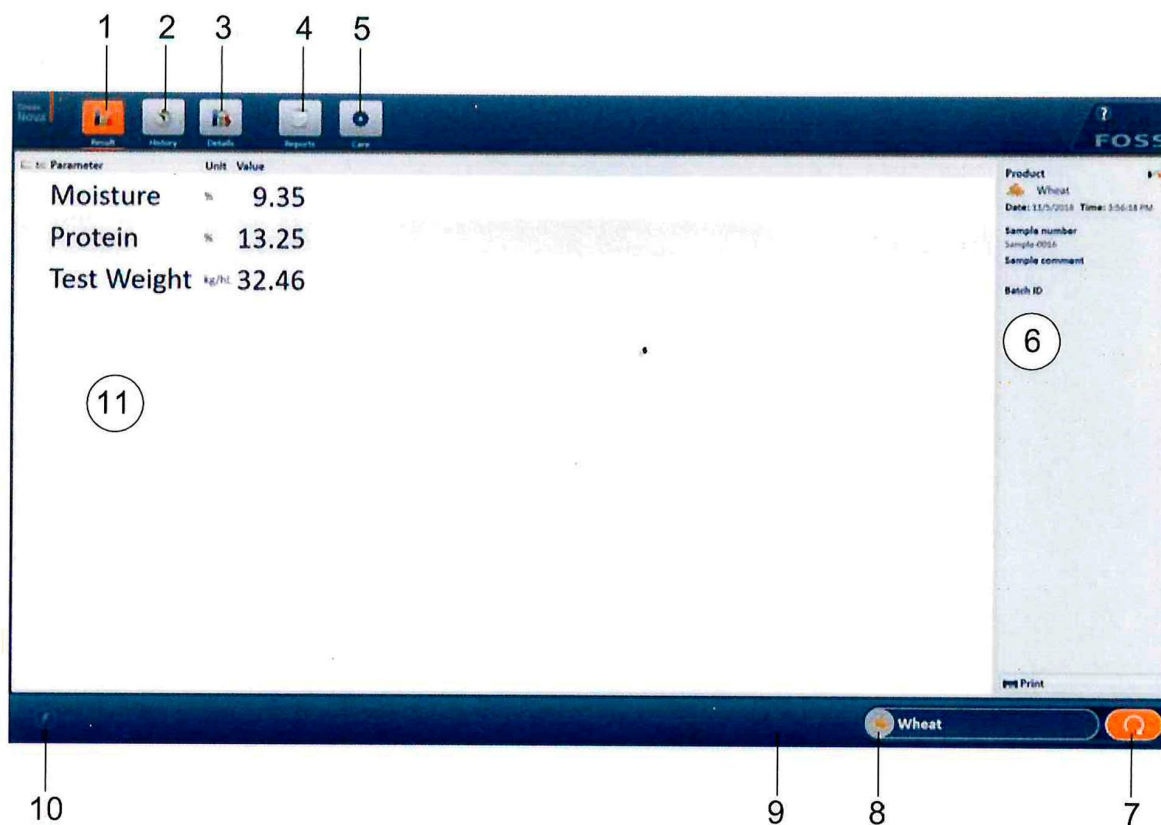


Fig. 1:1

- | | | | |
|---|---|----|---------------------|
| 1 | Results View | 7 | Start/Stop analysis |
| 2 | History View | 8 | Product Icon |
| 3 | Details | 9 | Progress |
| 4 | Reports | 10 | Event indicator |
| 5 | Care Menu | 11 | Result |
| 6 | Sample registration values
(User Defined Fields) | | |

Maintenance

See User Manual, chapter 5 for information about recommended maintenance.

Infratec™ - Daily Use


Start-Up

Switch on the instrument. Wait for the instrument start-up test to be completed.

Principle of Operation

The Infratec is a whole grain analyzer without need for sample preparation. Removal of stones, straws etc is recommended if the sample is contaminated with non-grain material. The path length is automatically adjusted according to the type of grain being analysed.

Analysing

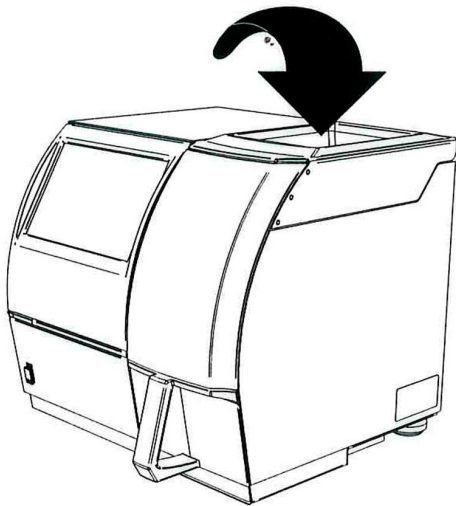1. Pour the sample in the hopper and press the start button on the screen.	
2. Select Product by pressing the Product Icon. Choose product to analyse from the Product list. Press Cancel if you do not want to change product.	